

Tobin Eyewash Bottles

An eyewash should be quick, simple and safe to use. The unique shape of Tobin's eyewash bottle insures that even a temporarily blinded and confused person can identify it as a safe eyewash.



Applications

Suitable when space is a problem or where several eyewash stations are required in the same area. Supplied with bottles, wall screws etc.

Features

- Provides 6 minutes washing time
- the bottle is made from blow-moulded plastic, filled with 1 litre of 0.9% saline, heat sealed, and then sterilised. It has a 3 year shelf life and remains sealed and sterile until use.
- the bottle hangs from a wall mounted rack by its cap. To open, simply break off the cap. An unopened bottle can be removed or replaced in the stand at will. An opened bottle cannot be replaced in the stand.
- the eye is washed by a soft flow consisting of six tiny streams. The gentle flow prevents chemicals or foreign bodies being driven further into damaged soft tissue
- the fact that the bottle is single use is a very important safety feature. Once the bottle has been opened, the liquid is no longer sterile. The round ended bottle is designed to prevent people saving unused liquid or refilling the bottle with other liquid. When empty the bottle has a hole at both ends, making it unsuitable for further use.
- tobins eyewash bottle has a round base and cannot stand alone. This is a particularly important safety feature that removes the temptation to take bottles from the stand and place them close to bottles with other contents. Mix-ups should never occur. It's important that everyone knows where the bottles are when an accident occurs.

TECHNICAL SPECIFICATIONS

Expiry date	Each bottle shows the recommended expiry date
Dimensions	390 x 200 x 100 mm



Tobin Eyewash Bottles
For a quick, simple and safe use